

13/11/19

Objectives

- learn to make NGM plates and pour them
- check on plates

Observations

- AML-238 SB - no embryos found on plate - could be sodium hydrochlorite at 14% is too strong - will be traditionally bleaching with Jun later in the week
- OG AML-283 ©
 - starved -> 15C
 - took about 6 days to reach this state - very very few crawled out of the chunk
 - picking practice
- (p) AML-283
 - also around 5-6 days until starved
 - picking practice
 - stored at 15C
- © AML-284 - ok - will probably be starved by Friday - can chunk for over the weekend
- (p) AML-284 x4 - numbers doubled in 48 hours or I picked more than 4 worms and now they're easier to see

From:
<http://wiki.scholz-lab.com/> - Scholz Lab Wiki

Permanent link:
http://wiki.scholz-lab.com/doku.php?id=wiki:documentation:wiki:documentation:nicolina:13_11_19&rev=1573659777

Last update: **2019/11/13 16:42**

