

Making a worm pick

1. Clip long glass end off to the point where it begins to fuse with the larger glass body
2. Cut to how long you want the wire to be - longer wires are easier for beginners (around 3-4cms)
3. Set up roaring bunsen
4. Put wire into glass -> heat both until glass wraps around the wire
5. Let cool
6. Flatten the tip of the wire as best as possible using a hard surface

From:

<http://wiki.scholz-lab.com/> - **Scholz Lab Wiki**

Permanent link:

http://wiki.scholz-lab.com/doku.php?id=wiki:documentation:wiki:documentation:nicolina:making_a_worm_pick

Last update: **2019/11/11 13:21**

